

# QIAseq Multimodal Panels

## Human Leukemia

**Cat. nos. 333932 UHS-009Z-12 and 333935 UHS-009Z-96**

**For consolidated targeted next-generation sequencing of DNA and RNA**


The QIAseq Multimodal Human Leukemia Panel permits the sequencing the whole coding region of 145 genes frequently mutated in leukemia as well as the detection of fusions to 176 genes also frequently observed in leukemia, with 6244 and 1116 primers respectively. Such frequently detected mutations and frequently detected fusions are likely to be relevant or even important to leukemia biology and may be used in translational research to classify samples with known phenotypes. Using the Catalogue of Somatic Mutations in Cancer (COSMIC), frequently mutated genes and fusions were collected for acute lymphoblastic leukemia (ALL), acute myeloid leukemia (AML), chronic lymphocytic leukemia (CLL), chronic myeloid leukemia (CML), chronic myelomonocytic leukemia (CMML), chronic neutrophilic leukemia (CNL), multiple myeloma, myelodysplastic syndromes (MDS), myeloproliferative neoplasms (MPN), other leukemias and other myeloid neoplasms. From the same cancer subtypes, additional fusions were also collected from the Archer Quiver Database, and more mutations and more fusions were both requested and suggested by QIAGEN KOLs and SMEs and found in other peer-reviewed publications. Using digital DNA sequencing, SNP/indels in, and fusions to, two different focused panels of genes relevant to leukemia can be detected with this primer pool.

For further details, consult the *QIAseq Multimodal Panels Handbook*.

### Shipping and storage

QIAseq Multimodal Panels (except QIAseq Beads, QIAseq Bead Binding Buffer, and Nuclease-free Water) are shipped on dry ice and should be stored immediately upon receipt at  $-30^{\circ}\text{C}$  to  $-15^{\circ}\text{C}$  in a constant-temperature freezer. QIAseq Beads, QIAseq Bead Binding Buffer, and Nuclease-free Water are shipped in a separate box on cold packs. Upon receipt, QIAseq Beads, QIAseq Bead Binding Buffer, and Nuclease-free Water should be stored at  $2-8^{\circ}\text{C}$ . QIAseq Multimodal Index kits are shipped on dry ice and should be stored at  $-30^{\circ}\text{C}$  to  $-15^{\circ}\text{C}$  upon arrival.

**Note:** Open the package and store the products appropriately immediately upon receipt.

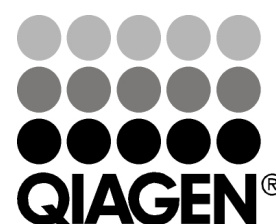

## Panel Contents

| QIAseq Multimodal Panel Catalog Number    | 333932                | 333935                 |
|-------------------------------------------|-----------------------|------------------------|
| Number of samples                         | 12                    | 96                     |
| Multimodal DHS Panel (DNA)                | 120 $\mu$ L           | 960 $\mu$ L            |
| Multimodal VHS Panel (RNA)                | 96 $\mu$ L            | 768 $\mu$ L            |
| Fragmentation Buffer, 10x                 | 40 $\mu$ L            | 192 $\mu$ L            |
| Fragmentation Enzyme Mix                  | 90 $\mu$ L            | 384 $\mu$ L            |
| FERA Solution                             | 15 $\mu$ L            | 60 $\mu$ L             |
| Side Reaction Reducer                     | 48 $\mu$ L            | 192 $\mu$ L            |
| FG Solution                               | 170 $\mu$ L           | 170 $\mu$ L            |
| DNA Ligase                                | 75 $\mu$ L            | 600 $\mu$ L            |
| Ligation Solution                         | 125 $\mu$ L           | 970 $\mu$ L            |
| Ligation Buffer, 5x                       | 160 $\mu$ L           | 1250 $\mu$ L           |
| DNA Ligation Adapter                      | 34 $\mu$ L            | 270 $\mu$ L            |
| ATP Solution                              | 36 $\mu$ L            | 290 $\mu$ L            |
| PAP Enzyme                                | 12 $\mu$ L            | 96 $\mu$ L             |
| PAP Dilution Buffer, 10x                  | 24 $\mu$ L            | 192 $\mu$ L            |
| T4 Polynucleotide Kinase                  | 12 $\mu$ L            | 125 $\mu$ L            |
| EZ Reverse Transcriptase                  | 15 $\mu$ L            | 150 $\mu$ L            |
| RNase Inhibitor                           | 12 $\mu$ L            | 96 $\mu$ L             |
| Multimodal RT Primer                      | 12 $\mu$ L            | 96 $\mu$ L             |
| Multimodal RT Buffer, 5x                  | 60 $\mu$ L            | 480 $\mu$ L            |
| Multimodal RT Enhancer                    | 12 $\mu$ L            | 48 $\mu$ L             |
| TEPCR Buffer, 5x                          | 60 $\mu$ L $\times$ 2 | 500 $\mu$ L $\times$ 2 |
| HotStarTaq® DNA Polymerase (6 U/ $\mu$ L) | 60 $\mu$ L            | 480 $\mu$ L            |
| Nuclease-free Water                       | 1 tube                | 10 mL                  |
| QIAseq Beads                              | 10 mL                 | 38.4 mL $\times$ 2     |
| QIAseq Bead Binding Buffer                | 10.2 mL               | 10.2 mL $\times$ 2     |

## Gene Table: DNA SNP/indel

| Gene Symbol | Gene Symbol | Gene Symbol | Gene Symbol |
|-------------|-------------|-------------|-------------|
| ABL1        | ADA         | ANKRD26     | ASXL1       |
| ASXL2       | ATM         | ATRX        | BCL6        |
| BCOR        | BCORL1      | BCR         | BIRC3       |
| BLM         | BRAF        | BRCA1       | BRCA2       |
| BRINP3      | C17orf97    | CALR        | CARD11      |
| CBL         | CBLB        | CBLC        | CDKN2A      |
| CEBPA       | CHEK2       | CREBBP      | CRLF2       |
| CSF1R       | CSF3R       | CTCF        | CUX1        |
| DAXX        | DDX41       | DNM2        | DNMT1       |
| DNMT3A      | EED         | EGFR        | ELANE       |
| EP300       | ETNK1       | ETV6        | EZH2        |
| FAM47A      | FAS         | FBXW7       | FLRT2       |
| FLT3        | GATA1       | GATA2       | GJB3        |
| GNA5        | HNRNPBK     | HRA5        | IDH1        |
| IDH2        | IKZF1       | IKZF3       | IL7R        |
| JAK1        | JAK2        | JAK3        | KAT6A       |
| KCNA4       | KCNK13      | KDM6A       | KDR         |
| KIT         | KLHDC8B     | KLHL6       | KMT2A       |
| KMT2C       | KMT2D       | KRAS        | LRRC4       |
| LUC7L2      | MAP2K1      | MLH1        | MPL         |
| MSH2        | MSH6        | MYC         | MYD88       |
| NBN         | NF1         | NOTCH1      | NPAT        |
| NPM1        | NRAS        | NSD1        | NTRK3       |
| OR13H1      | OR8B12      | P2RY2       | PAX5        |
| PBRM1       | PCDHB1      | PDGFRA      | PHF6        |
| PML         | PMS2        | PPM1D       | PRAMEF2     |
| PRF1        | PRPF40B     | PRPF8       | PTEN        |
| PTPN11      | RAD21       | RB1         | RELN        |
| RET         | RUNX1       | SAXO2       | SETBP1      |
| SF1         | SF3A1       | SF3B1       | SH2B3       |
| SH2D1A      | SMARCB1     | SMC1A       | SMC3        |
| SRP72       | SRSF2       | STAG2       | STAT3       |
| STXBP2      | SUZ12       | TAL1        | TERC        |
| TERT        | TET2        | TNFRSF13B   | TP53        |
| TPMT        | TUBA3C      | U2AF1       | U2AF2       |
| WAS         | WRN         | WT1         | XPO1        |
| ZRSR2       |             |             |             |

## Gene Table: RNA Fusions

| Gene Symbol | Gene Symbol | Gene Symbol | Gene Symbol |
|-------------|-------------|-------------|-------------|
| ABLI        | ABLI2       | ABLI        | ABLI2       |
| ACSL6       | ACTN4       | AFDN        | AFF1        |
| AFF3        | AFF4        | AGGF1       | ALK         |
| ANLN        | ARHGAP26    | ARHGEF12    | ASNS        |
| ATF7IP      | AUTS2       | BCL9        | BCOR        |
| BCR         | BTBD18      | CASC5       | CBFA2T3     |
| CBFB        | CBL         | CCDC88C     | CEP170B     |
| CHD1        | CHIC2       | CREBBP      | CRLF2       |
| CSF1R       | CT45A2      | DAB2IP      | DDX10       |
| DDX3X       | DEK         | DHH         | EBF1        |
| ELL         | ELN         | EML1        | EP300       |
| EP515       | ERG         | ESRRB       | ETV6        |
| EWSR1       | FGFR1       | FGFR1OP     | FIP1L1      |
| FLNA        | FLT3        | FNBP1       | FOXO3       |
| FOXO4       | FOXP1       | FRYL        | FUS         |
| GAS7        | GATA1       | GLIS2       | GPHN        |
| HAUS1       | HLF         | HNRNPUL1    | HOXA9       |
| HOXC11      | HOXC13      | HOXD13      | IKZF3       |
| INPP5D      | IRF2BP2     | JAK2        | KAT6A       |
| KDM5A       | KIAA1524    | KMT2A       | KMT2E       |
| LASP1       | LNP1        | LPP         | LYN         |
| MAML2       | MAPRE1      | MECOM       | MEF2D       |
| MKL1        | MLF1        | MLLT1       | MLLT10      |
| MLLT11      | MLLT3       | MLLT6       | MN1         |
| MNX1        | MYB         | MYC         | MYH11       |
| MYO1F       | NCKIPSD     | NCOA2       | NDE1        |
| NFI         | NOTCH1      | NPM1        | NSD1        |
| NTRK3       | NUMA1       | NUP214      | NUP98       |
| P2RY8       | PAX5        | PBX1        | PCM1        |
| PDCD1LG2    | PDGFRA      | PDGFRB      | PDS5A       |
| PHF23       | PICALM      | PML         | PRDM16      |
| PRKG2       | PRRX1       | PSIP1       | PTPRR       |
| RANBP2      | RAP1GDS1    | RARA        | RBM15       |
| RCS1        | RHEBL1      | ROS1        | RPL22       |
| RPN1        | RUNX1       | RUNX1T1     | SALL2       |
| SARNP       | SEPT2       | SEPT5       | SEPT6       |
| SEPT9       | SET         | SETD2       | SFPQ        |
| SH3GL1      | SORBS2      | SPI1        | SSBP2       |
| STAT5B      | STIL        | STMN1       | TAF15       |
| TAL1        | TCF3        | TCF7        | TET1        |
| TFG         | TNIP1       | TOP1        | TOP3A       |
| TRIP11      | UBA2        | USP42       | USP9X       |
| WH5C1L1     | WTIP        | ZBTB16      | ZFYVE19     |
| ZMIZ1       | ZMYM2       | ZNF384      | ZNF521      |

## Functional Groupings

### DNA SNP/INDELS:

**Acute Lymphoblastic Leukemia (ALL):** ASXL2, ATM, BRAF, CALR, CDKN2A, CREBBP, CRLF2, CSF3R, CTCF, DNM2, EGFR, EP300, FBXW7, GATA2, HNRNPK, HRAS, IKZF3, IL7R, KDM6A, KDR, KMT2C, LRRC4, MAP2K1, MLH1, MSH2, MSH6, NOTCH1, NTRK3, PAX5, PDGFRA, PMS2, PRAMEF2, PTEN, RELN, SMARCB1.

**Acute Myeloid Leukemia (AML):** ANKRD26, ASXL1, ATM, BCOR, BCORL1, BIRC3, BRAF, C17orf97, CALR, CARD11, CBLC, CDKN2A, CEBPA, CHEK2, CREBBP, CSF1R, CSF3R, CTCF, DAXX, DDX41, DNM2, DNMT1, ELANE, EP300, FLRT2, FLT3, GATA1, GATA2, HNRNPK, IDH1, IDH2, IKZF1, IL7R, JAK1, JAK3, KDM6A, KDR, KIT, KMT2A, KMT2C, KMT2D, KRAS, LRRC4, MAP2K1, MPL, MSH6, MYC, NBN, NOTCH1, NPM1, NRAS, NSD1, NTRK3, OR13H1, OR8B12, P2RY2, PBRM1, PCDHB1, PDGFRA, PHF6, PPM1D, PRAMEF2, PRPF8, PTEN, PTPN11, RAD21, RET, RUNX1, SF1, SF3A1, SMARCB1, SMC1A, SMC3, SRP72, SRSF2, STAG2, STXBP2, U2AF1, U2AF2, WT1.

**Chronic Lymphocytic Leukemia (CLL):** ADA, BIRC3, BLM, BRAF, CALR, CHEK2, CSF3R, KCNA4, KLHL6, KMT2C, MAP2K1, NBN, NPAT, NTRK3, OR13H1, OR8B12, PRAMEF2, SRP72, TAL1, TERT, TUBA3C, WAS, WRN.

**Chronic Myeloid Leukemia (CML):** ABL1, CALR, CDKN2A, CEBPA, CREBBP, CSF1R, CSF3R, FBXW7, GATA2, KDM6A, MSH2, MSH6, RB1, SMC1A, TP53.

**Chronic Myelomonocytic Leukemia (CMML):** CALR, CEBPA, CSF1R, CSF3R, HRAS, KMT2C, LUC7L2, SRSF2.

**Chronic Neutrophilic Leukemia (CNL):** CALR, CSF3R.

**Multiple Myeloma:** ATM, BCL6, BCR, BIRC3, BRAF, CDKN2A, CEBPA, EGFR, FBXW7, GJB3, HRAS, KDM6A, MYC, NOTCH1, PTEN, SH2D1A, SMARCB1.

**Myelodysplastic Syndromes (MDS):** ATRX, CALR, CDKN2A, CEBPA, CSF1R, CSF3R, EP300, ETNK1, GNAS, HRAS, KDM6A, KMT2A, KMT2C, RAD21, RB1, SETBP1, SF1, SF3A1, SMC3, SRSF2, STAG2, U2AF1, U2AF2, XPO1, ZRSR2.

**Myeloid Malignancies:** CBL, CBLB, DNMT3A, EED, ETV6, EZH2, PRPF40B, SUZ12, TET2, TP53.

**Myeloproliferative Neoplasms (MPN):** ABL1, ASXL1, CALR, CSF1R, JAK2, JAK3, KAT6A, KRAS, MPL, NF1, NRAS, RB1, SETBP1, SF3B1, SH2B3, SRSF2, STAG2.

**Myelofibrosis:** CALR, CHEK2, CSF1R, IDH1, IDH2, SRSF2.

**Other Myeloid Neoplasms:** BRAF, CDKN2A, CEBPA, FBXW7, HRAS, IKZF3, KLHDC8B, KMT2C, MSH6, NTRK3, PTEN, SRP72, TPMT.

**Other Myeloid Neoplasm Genes:** BRCA1, BRCA2, BRINP3, CUX1, FAM47A, FAS, KCNK13, MYD88, PML, PRF1, SAXO2, STAT3, TERC, TNFRSF13B.

---

## **GENE FUSIONS:**

**Acute Leukemia of Ambiguous Lineage:** ABL1, AFF1, BCR, CT45A2, ETV6, EWSR1, GPHN, KMT2A, MLLT10, MN1, NUP214, PBX1, SET, TCF3, ZNF384.

**Acute Lymphoblastic B Cell Leukemia:** ABL1, ACTN4, AFF1, AFF3, AFF4, BCR, BTBD18, EPS15, ERG, ETV6, EWSR1, FUS, GAS7, JAK2, KMT2A, MAPRE1, MLLT1, NUP214, PAX5, PBX1, PCM1, RUNX1, SET, SSBP2, TCF3, ZNF384.

**Acute Lymphoblastic Leukemia (ALL):** ABL1, ABL2, ACTN4, AFF1, AFF3, AFF4, AGGF1, ASNS, ATF7IP, AUTS2, BCL9, BCR, BTBD18, CASC5, CREBBP, CRLF2, CSF1R, DDX3X, EBF1, ELL, ELN, EML1, EP300, EPS15, ESRRB, ETV6, EWSR1, FGFR1, FGFR1OP, FLT3, FOXO4, FOXP1, FRYL, GAS7, GPHN, HLF, HNRNPUL1, INPP5D, JAK2, KMT2A, KMT2E, LNP1, MAML2, MEF2D, MLLT1, MLLT10, MLLT3, AFDN, MN1, MNX1, MYC, NCOA2, NOTCH1, NUP214, NUP98, P2RY8, PAX5, PBX1, PCM1, PDGFRB, PICALM, PML, PRDM16, RAPIGDS1, RARA, RCSD1, RUNX1, SALL2, SET, SFPQ, SPI1, SSBP2, STIL, STMN1, TAF15, TAL1, TCF3, TCF7, TET1, USP9X, ZBTB16, ZMIZ1, ZMYM2, ZNF384, ZNF521.

**Acute Lymphoblastic T Cell Leukemia:** ABL1, BCR, CASC5, ETV6, JAK2, KMT2A, MLLT1, MLLT10, AFDN, NOTCH1, NUP214, PBX1, RUNX1, SET, STIL, TAL1, TCF3.

**Acute Monoblastic Leukemia:** KMT2A, MLLT10, PICALM.

**Acute Myeloid Leukemia (AML):** ABI1, ABI2, ABL1, AFF1, AFF4, ALK, ANLN, ARHGAP26, ARHGEF12, BCOR, BCR, CASC5, CBFA2T3, CBL, CCDC88C, CEP170B, CHD1, CHIC2, CREBBP, DAB2IP, DDX10, DEK, DHH, ELL, EP300, EPS15, ERG, ETV6, FGFR1, FIP1L1, FLNA, FNBP1, FOXO3, FOXO4, FRYL, FUS, GAS7, GATA1, GLIS2, GPHN, HAUS1, HOXA9, HOXC11, HOXC13, HOXD13, IRF2BP2, JAK2, KAT6A, KDM5A, KIAA1524, KMT2A, LASP1, LNP1, LPP, LYN, MAML2, MECOM, MKL1, MLF1, MLLT1, MLLT10, MLLT11, MLLT3, AFDN, MLLT6, MN1, MNX1, MYB, MYH11, MYO1F, NCKIP5D, NCOA2, NF1, NPM1, NSD1, NTRK3, NUMA1, NUP214, NUP98, PBX1, PCM1, PDGFRA, PDGFRB, PDS5A, PHF23, PICALM, PML, PRDM16, PRRX1, PSIP1, PTPRR, RANBP2, RARA, RBM15, RHEBL1, RPL22, RPN1, RUNX1, RUNX1T1, SARNP, SEPT2, SEPT5, SEPT6, SEPT9, SET, SH3GL1, SORBS2, STAT5B, TAF15, TCF3, TET1, TFG, TOP1, TOP3A, TRIP11, UBA2, USP42, WHSC1L1, WTIP, ZBTB16, ZFYVE19, ZNF384.

**B Cell Prolymphocytic Leukemia:** MYC.

**Burkitt Leukemia:** MYC.

**Chronic Lymphocytic Leukemia (CLL):** MYC, NOTCH1.

**Chronic Myeloid Leukemia (CML):** ABL1, BCR, CCDC88C, ETV6, JAK2, MECOM, MN1, NDE1, PCM1, PDGFRB, RUNX1.

**Chronic Myelomonocytic Leukemia (CMML):** CREBBP, ELL, ETV6, KMT2A, MN1.

**Lymphoblastic Leukemia:** AFF1, KMT2A.

**Myelodysplastic Syndromes (MDS):** ABL1, ACTN4, ALK, CASC5, CREBBP, DEK, ERG, ETV6, FRYL, FUS, HOXD13, JAK2, KMT2A, LNP1, MAML2, MECOM, MLF1, MN1, MNX1, NDE1, NPM1, NUP214, NUP98, PCM1, PDGFRB, PRDM16, PTPRR, RANBP2, RUNX1, RUNX1T1, TOP1.

**Myeloproliferative Neoplasms (MPN):** ABL1, ACSL6, BCR, ETV6, FGFR1, FGFR1OP, FIP1L1, FLT3, HOXA9, HOXD13, JAK2, LYN, MECOM, MLF1, MN1, NPM1, NUP98, PDGFRB, PSIP1, RARA, RUNX1, TNIP1, ZMYM2.

**Plasma Cell Leukemia:** MYC.

**T Cell Prolymphocytic Leukemia:** MYC.

**T-Cell Acute Myelocytic Leukemia:** KMT2A, MLLT10.

**Other Leukemias:** IKZF3, PDCD1LG2, PRKG2, ROS1, SETD2.

---

## Ordering Information

| Product                             | Contents                                                                                                                                                                                                                                                                  | Cat. no. |
|-------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|----------|
| QIAseq Multimodal Panel (12)        | Kit containing all reagents (except indexes) for multimodal (DNA and RNA) sequencing; fixed small panel for a total of 12 samples: 12 DNA and 12 RNA libraries                                                                                                            | 333932   |
| QIAseq Multimodal Panel (96)        | Kit containing all reagents (except indexes) for multimodal (DNA and RNA)                                                                                                                                                                                                 | 333935   |
| QIAseq Multimodal HC Panel (12)     | Kit containing all reagents (except indexes) for multimodal (DNA and RNA) sequencing; fixed high content (HC) panel for a total of 12 samples: 12 DNA and 12 RNA libraries                                                                                                | 333942   |
| QIAseq Multimodal HC Panel (96)     | Kit containing all reagents (except indexes) for multimodal (DNA and RNA) sequencing; fixed high content (HC) panel for a total of 96 samples: 96 DNA and 96 RNA libraries                                                                                                | 333945   |
| QIAseq Multimodal Custom Panel (96) | Kit containing all reagents (except indexes) for multimodal (DNA and RNA) sequencing; custom panel for a total of 96 samples: 96 DNA and 96 RNA libraries                                                                                                                 | 333955   |
| QIAseq Multimodal Index I SW (12)   | Box containing oligos, enough to process a total of 12 samples, for indexing up to a                                                                                                                                                                                      | 333982   |
| QIAseq Multimodal HT SW 96 UDI      | Box containing oligos, enough to process a total of 96 samples, for indexing up to a total of 96 samples (96 for DNA and 96 for RNA) for multimodal panel sequencing on Illumina platforms. Note: For Separated Enrichment Workflow (SW) for Targeted Enrichment          | 333986   |
| QIAseq Multimodal Index I (12)      | Box containing oligos, enough to process a total of 12 samples, for indexing up to a total of 12 samples (12 for DNA and 12 for RNA libraries) for multimodal panel sequencing on Illumina platforms. Note: For Combined Enrichment Workflow (CW) for Targeted Enrichment | 333962   |
| QIAseq Multimodal HT CW 96 UDI      | Box containing oligos, enough to process a total of 96 samples, for indexing up to a                                                                                                                                                                                      | 333979   |
| <b>Related products</b>             |                                                                                                                                                                                                                                                                           |          |
| QIAseq DNA QuantiMIZE Assay Kit     | qPCR assays for optimizing amount of input DNA and PCR cycling conditions for targeted enrichment of FFPE DNA                                                                                                                                                             | 333414   |
| QIAamp® DNA Mini Kit (50)           | For 50 DNA preps: 50 QIAamp Mini Spin Columns, QIAGEN Proteinase K, collection tubes (2 mL), reagents and buffers                                                                                                                                                         | 51304    |

| Product                                  | Contents                                                                                                                                                        | Cat. no. |
|------------------------------------------|-----------------------------------------------------------------------------------------------------------------------------------------------------------------|----------|
| QIAamp Circulating Nucleic Acid Kit (50) | For 50 DNA preps: QIAamp Mini Columns, Tube Extenders (20 mL), QIAGEN Proteinase K, carrier RNA, buffers, VacConnectors, and collection tubes (1.5 mL and 2 mL) | 55114    |
| AllPrep DNA/RNA Mini Kit (50)            | For 50 minipreps: AllPrep DNA Spin Columns, RNeasy Mini Spin Columns, collection tubes, RNase-free water, and buffers                                           | 80204    |
| AllPrep DNA/RNA FFPE Kit (50)            | 50 RNeasy MinElute Spin Columns, 50 QIAamp MinElute Spin Columns, collection tubes, RNase-free reagents, and buffers                                            | 80234    |
| PAXgene Blood DNA Kit (25)               | Processing tubes and buffers for 25 preparations                                                                                                                | 761133   |
| PAXgene Blood RNA Kit (50)               | 50 PAXgene Spin Columns, 50 PAXgene Shredder Spin Columns, processing tubes, RNase-Free DNase I, RNase-free reagents and buffers                                | Inquire  |

QIAseq Multimodal Panels are intended for molecular biology applications. These products are not intended for the diagnosis, prevention or treatment of a disease.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at [www.qiagen.com](http://www.qiagen.com) or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN® (QIAGEN Group)

1100784 02/2016 © 2016 QIAGEN, all rights reserved.

[www.qiagen.com](http://www.qiagen.com)

Australia ■ 1-800-243-800

Austria ■ 0800/281010

Belgium ■ 0800-79612

Brazil ■ 0800-557779

Canada ■ 800-572-9613

China ■ 8621-3865-3865

Denmark ■ 80-885945

Finland ■ 0800-914416

France ■ 01-60-920-930

Germany ■ 02103-29-12000

Hong Kong ■ 800 933 965

Ireland ■ 1800 555 049

Italy ■ 800-787980

Japan ■ 03-6890-7300

Korea (South) ■ 080-000-7145

Luxembourg ■ 8002 2076

Mexico ■ 01-800-7742-436

The Netherlands ■ 0800 0229592

Norway ■ 800-18859

Singapore ■ 1800-742-4368

Spain ■ 91-630-7050

Sweden ■ 020-790282

Switzerland ■ 055-254-22-11

UK ■ 01293-422-911

USA ■ 800-426-8157



## Sample to Insight